

## SEPARATION OF METHYLCLOTHIAZIDE USING NEW IMMOBILIZED COLUMN – CHIRALPAK IG

### APPLICATION NOTE

#### INTRODUCTION

Daicel Corporation has introduced a new and unique chiral selector, CHIRALPAK® IG, to its line of Daicel immobilized chiral stationary phases (CSPs). The IG phase is an amylose polymer derivatized with phenyl moieties (chloro- and methyl-groups) in meta positions. This distinctive phase was developed from our extensive library of > 150 CSPs. Moreover, our studies have shown that the IG chiral selector has the highest selectivity of any of our immobilized CSPs.

Meta-substituted immobilized chiral selectors such as CHIRALPAK IA, IB and IC have been shown to have remarkable affinity for resolution of chiral compounds from different types of molecules. Based on laboratory results, we have found that the addition of CHIRALPAK IG forms the most effective set of four chiral columns, CHIRALPAK IA, IB, IC, and IG, to be utilized for primary screening. It is noteworthy that the four best columns are all meta-substituted.

#### Primary Screening Columns

- CHIRALPAK IA: Amylose tris (3,5-dimethylphenylcarbamate)
- CHIRALPAK IB: Cellulose tris (3,5-dimethylphenylcarbamate)
- CHIRALPAK IC: Cellulose tris (3,5-dichlorophenylcarbamate)
- CHIRALPAK IG: Amylose tris (3-chloro-5-methylphenylcarbamate)

#### EXPERIMENTAL AND DISCUSSION

A CHIRALPAK IG column, packed with 5- $\mu$ m particles was used to develop the enantioselective separation of methyclothiazide. The mobile phase was a mixture of ammonium bicarbonate and acetonitrile. All chromatographic conditions are given in Figure 1.

#### SEPARATION EXAMPLE – REVERSED PHASE

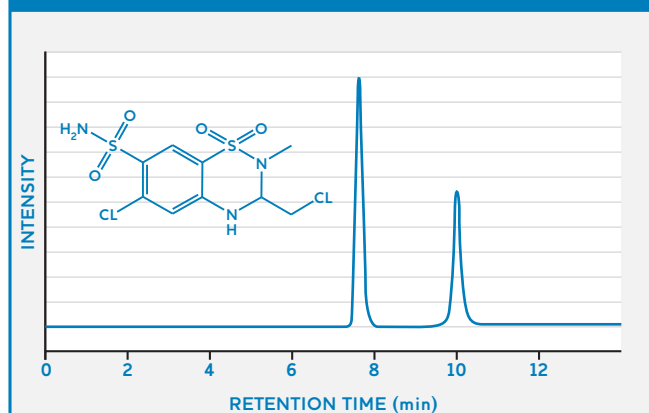


Figure 1: Separation of Methyclothiazide

#### CHROMATOGRAPHIC CONDITIONS

<b>Column Size:</b>	Daicel CHIRALPAK IG 4.6 mm i.d. x 250 mm long
<b>Mobile Phase:</b>	20mM $\text{NH}_4\text{HCO}_3$ aq/MeCN (pH 9 with DEA, 60/40)
<b>Flow Rate:</b>	1 ml/min
<b>UV Detection:</b>	254 nm
<b>Column Temperature:</b>	25 °C

**Note:** Methyclothiazide is a thiazide diuretic that is used to treat high blood pressure (hypertension). Using CHIRALPAK IG provides fast and effective separation of methyclothiazide enantiomers



#### CHIRAL TECHNOLOGIES, INC.

800 North Five Points Road  
West Chester, PA 19380 USA  
Tel: 610-594-2100  
Fax: 610-594-2325  
www.chiraltech.com  
email: chiral@chiraltech.com

#### CHIRAL TECHNOLOGIES EUROPE

Parc d'Innovation, Bd Gonthier d'Andernach  
67404 Illkirch Cedex, France  
Tel: +33 (0) 388 79 52 00  
Fax: +33 (0) 388 66 71 66  
www.chiraltech.com  
e-mail: cte@chiral.fr

