

CHIRAL TECHNOLOGIES INC

DAICEL GROUP

Innovation Begins Here

On the method development of immobilized polysaccharide chiral stationary phases in SFC using extended range of co-solvents

William L. Watts¹, Jr., J.T. Lee¹, James Lee¹, Joseph Barendt¹, Frank Riley² and Pilar Franco³

- ¹ CHIRAL TECHNOLOGIES INC.
- ² CHIRAL TECHNOLOGIES EUROPE
- ³ PFIZER

SFC 2013



SFC as efficient screening tool for chiral molecules

PRIMARY SCREENING

Limited selection of Chiral Stationary Phases (CSPs)


Co-solvents

EtOH, MeOH, 2-PrOH

(sometimes Acetonitrile)

Approach with high success rate!

Broadly accepted

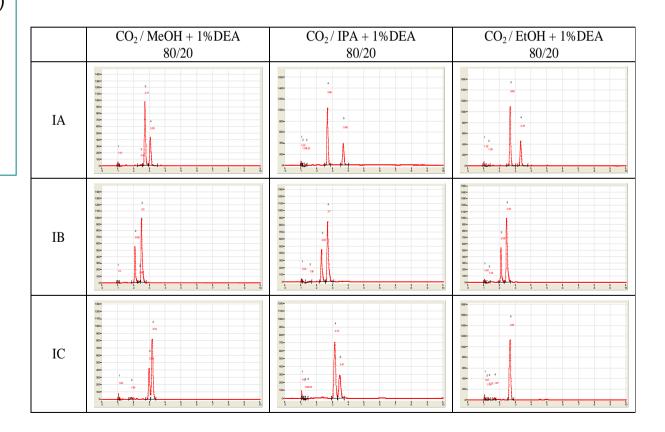



How choosing the CSPs?

Broad recognition

Most general purpose screening

Analytical or preparative application

Limited number of columns




Polysaccharide-derived selectors




A number of CSPs available

RO	Amylose-based		
	CSP	Nature	-R
	CHIRALPAK AD-H	Coated	H—CH ₃

CHIRALPAK IA	Immobilized	O CH ₃
CHIRALPAK ID	Immobilized	N CI
CHIRALPAK IE	Immobilized	O CI
CHIRALPAK IF	Immobilized	$ N$ $ CH_3$

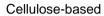

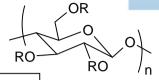

CSP	Nature	-R
CHIRALCEL OD-H	Coated	O CH ₃
CHIRALCEL OJ-H	Coated	O CH_3
CHIRALCEL OZ-H	Coated	O CI
CHIRALCEL OX-H	Coated	$ \begin{array}{c} H \\ O \\ CH_3 \end{array} $

CHIRALPAK IB	Immobilized	$\begin{array}{c} H \\ N \\ O \\ CH_3 \end{array}$
CHIRALPAK IC	Immobilized	H N O CI

Screening works but there are a number of aspects to be investigated



The start of the project

- ➤ CTI and Pfizer Pharmaceuticals in Groton started the first discussions about the project in 2010
- > A joint project was setup to learn more about:
 - > Screening strategies for the resolution of enantiomers in SFC
 - > The use of extended solvent range
 - > Additional options offered for prep use
- Pfizer loaned an analytical SFC to CTI for the project



Why extended solvent range in SFC could be of interest?

- ✓ Solubility reasons
- ✓ Stability issues
- ✓ Increase of screening success rate



Sample solubility




CHIRALPAK IC (250 x 30 mm) CO₂/EtOH 70/30 120 ml/min, 25℃


DCM is less polar than alcohols or THF (not compatible with all columns)

Solubility in EtOH < 2 g/L Solubility in EtOH/DCM 90/10 = 58 g/L

Injecting in a solvent different from mobile phase

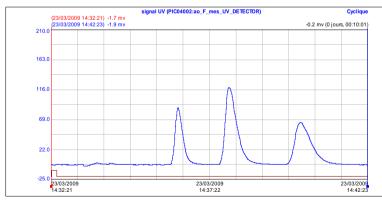

Analytical injection

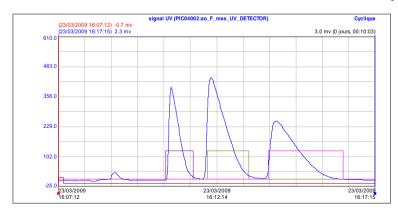



Injection in EtOH/DCM 90/10 - 2ml - 116 mg



Sample stability

A case study in LC

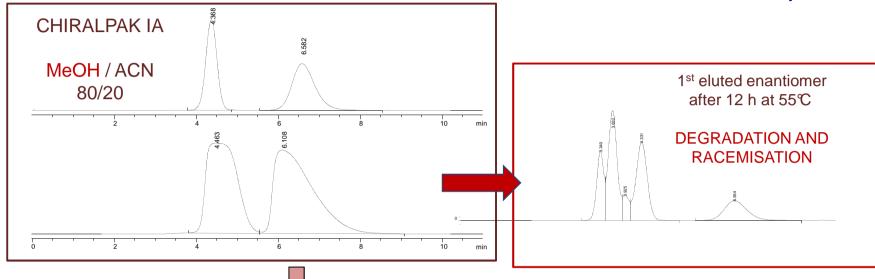

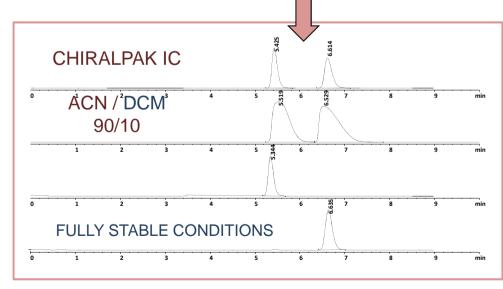

Compound with glutarimide moiety

In certain cases, degradation and/or racemisation can be controlled with lower evaporation temperature




Extended Range Co-Solvent System

- ✓ Evaluations of ethyl acetate, THF and DCM
- ✓ Use of immobilized-type polysaccharide CSPs
- ✓ Parameters investigated:
 - ✓ Elution strength of those solvents in SFC
 - ✓ Role of the addition of alcohol in the elution and the success rate
 - ✓ Role of additives for the elution and resolution of acidic and basic molecules



Structure of CSPs used in the study

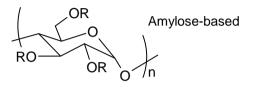

CSP	Nature	-R
CHIRALPAK IA	Immobilized	$ \begin{array}{c} $
CHIRALPAK ID	Immobilized	H N O Cl
CHIRALPAK IE	Immobilized	H—————————————————————————————————————
CHIRALPAK IF*	Immobilized	O CI

-R	
CH_3 CH_3	
CI	

^{*} Not available when study was performed

Gradient or isocratic?

Journal of Chromatography B, 875 (2008) 230-236



Contents lists available at ScienceDirect

Journal of Chromatography B

journal homepage: www.elsevier.com/locate/chromb



Preparative chromatographic resolution of racemates using HPLC and SFC in a pharmaceutical discovery environment[☆]

Larry Miller*, Matt Potter

Amgen, One Kendall Square, Bldg 1000, Cambridge, MA 02139, USA

ARTICLE INFO

Article history: Received 18 April 2008 Accepted 17 June 2008 Available online 1 July 2008

Keywords: Chiral separations Preparative chromatography Supercritical fluid chromatography Chiral stationary phase

ABSTRACT

The preparative chromatographic resolution of racemates has become a standard approach for the generation of enantiomers in pharmaceutical discovery laboratories. This paper will discuss the use of preparative HPLC and SFC to generate individual enantiomers for discovery activities. Analytical HPLC and SFC method development to rapidly screen chiral stationary phases and solvent combinations will be presented. The usefulness of preparative chromatographic resolution of racemates will be demonstrated through the presentation of numerous non-routine case studies from the laboratories at Amgen.

© 2008 Elsevier B.V. All rights reserved.

L. Miller et al., J. Chromatogr. B 875 (2008) 230-236

- Combination of LC and SFC
- Initial screening in gradient mode
- 3. Final isocratic method

Literature example





Gradient or isocratic?

Some examples in SFC


Back pressure: 150bar, Flow rate: 3.0ml/min,

Temperature: 35°C

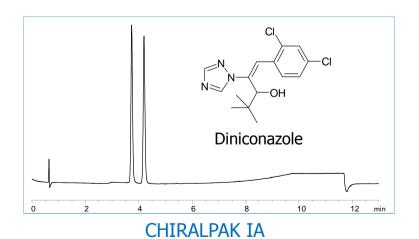

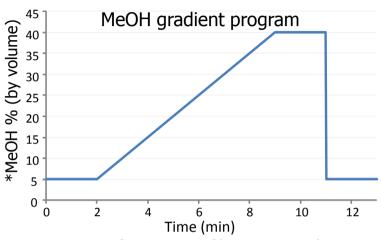


* Containing 0.3% DEA for screening of basic compounds

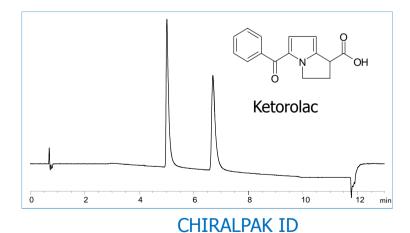



From gradient to isocratic

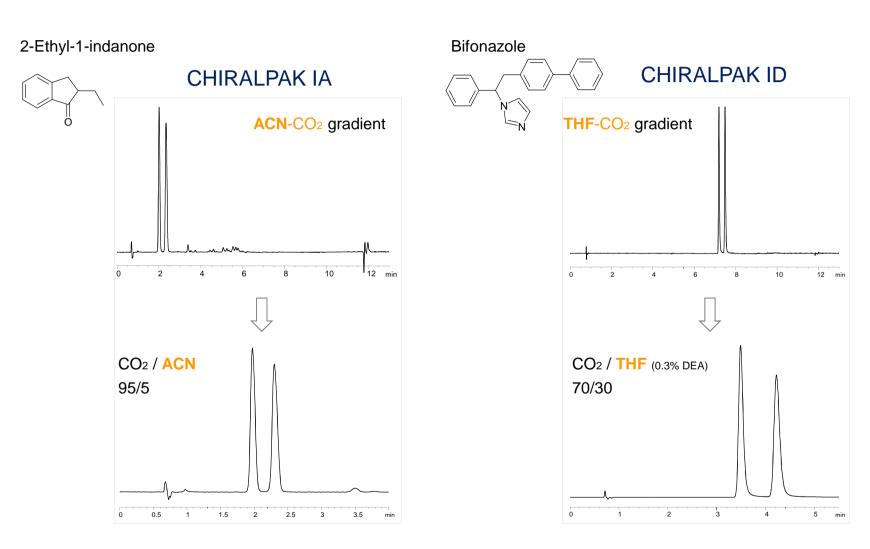

Easy transfer of conditions



SFC Conditions used in the study

System	Berger Analytical System
Column	IA, IB, IC, ID, IE
Column Size	250 x 4.6 mm, 5 μm
Mobile	A: CO ₂ , B: Modifier (co-solvents)
Phases	See next slide
Gradient	5 – 50% B gradient at 6.5%/min, hold for 3 mins, re-equilibrate for 2 mins
Temp.	40° C
Pressure	120 bar
Detection	DAD: 215, 230, 254 nm
Flow	4.0 mL/min
Injection	10 μL
Run Time	12 min



Evolution screening conditions in the study

Initial setup

- 100% THF
- 100% EtOAc
- 50% DCM/50% MeOH
- 0.1% IPAmine added to co-solvents for basic molecules
- No acid additive for acidic molecules

Revised setup

- 95% THF/5% MeOH
- 90% EtOAc/10% MeOH
- 90% DCM/10% MeOH



• 0.2% IPAmine added to co-solvents for basic molecules



• 0.2% TFA added to cosolvents for acidic molecules



Selected Acidic Molecules

Code #	ACIDS
acid 1	trans-2-Phenyl-1-cyclopropane carboxylic acid
acid 2	Bis-N-Cbz-DL-Lysine
acid 3	Sulindac
acid 4	Suprofen
acid 5	Proglumide
acid 6	Dichlorprop
acid 7	Fenoprofen, Ca-Salt
acid 8	2-(4-Hydroxy-phenyl)-propionic acid
acid 9	3-Phenylbutyric acid
acid 10	2-Phenylbutyric acid
acid 11	2-Phenylpropionic acid

11 acids



Selected Basic Molecules


Sample Code	BASES
basic 1	Mebeverine HCl
basic 2	Propafenone HCl
basic 3	Homatropine
basic 4	Labetalol HCl
basic 5	Verapamil HCl
basic 6	Chlophedianol HCl
basic 7	Acebutolol
basic 8	Atropine
basic 9	Ofloxacin (zwitterionic)
basic 10	Diperodon HCl
basic 11	Ketamine HCI
basic 12	Nadolol
basic 13	Salbutamol HCl
basic 14	p-Chloroamphetamine HCl

14 bases




The elution strength

THF mixtures

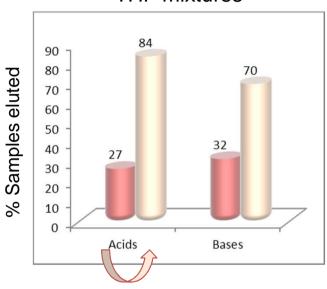


Addition of 10% MeOH to THF

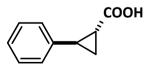

Trans-2-phenyl-1-cyclopropane carboxylic acid

CHIRALPAK IA

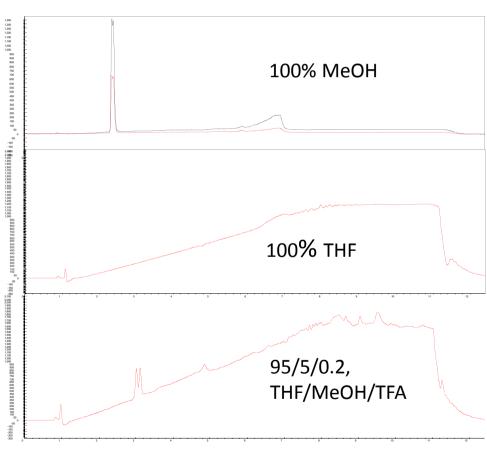




The elution strength


Percentages of compounds eluted

THF mixtures

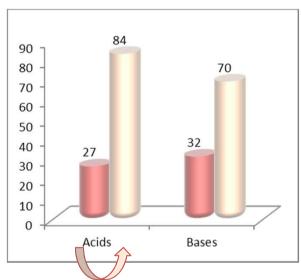


Addition of 10% MeOH to THF

EtOAc mixtures

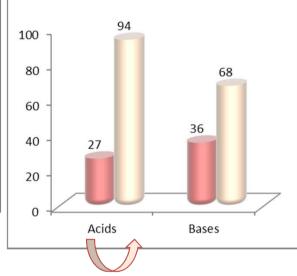

Addition of 10% MeOH to EtOAc

DCM mixtures

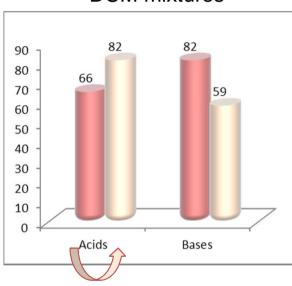

From 50% MeOH to 10% MeOH in DCM

Statistics on all runs and conditions tested (total: 125)




Bringing selectivity with extended solvent range


	ACIDS			
		Alcohols	Old Ext. run	New Ext. run
acid 1	trans-2-Phenyl-1-cyclopropane carboxylic acid	8	4	10
acid 2	Bis-N-Cbz-DL-Lysine	6	none eluted	3
acid 3	Sulindac	6	2	4
acid 4	Suprofen	8	4	12
acid 5	Proglumide	11	4	12
acid 6	Dichlorprop	2	3	8
acid 7	Fenoprofen, Ca-Salt	0 resolved	3	5
acid 8	2(4-HO-PheO)-propionic acid	5	4	13
acid 9	3-Phenylbutyric acid	1	1	6
acid 10	2-Phenylbutyric acid	0 resolved	0 resolved	4
acid 11	2-Phenylpropoionic acid	0 resolved	0 resolved	3
	BASES			
		Alcohols	Old Ext. run	New Ext. run
basic 1	Mebeverine HCI	9	4	6
basic 2	Propafenone HCI	9	5	7
basic 3	Homatropine	7	1	2
basic 4	Labetalol HCI	5	1	4
basic 5	Verapamil HCI	4	3	7
basic 6	Chlophedianol HCI	6	5	10
basic 7	Acebutolol	4	1	2
basic 8	Atropine	6	3	1 (IC)
basic 9	Ofloxacin	2	none eluted	1 (ID)
basic 10	Diperodon HCI	11	10	13
basic 11	Ketamin HCI	2	5	8
basic 12	Nadolol	12	3	1 (IA)
basic 13	Salmbutamol HCI	4	0 resolved	1 (IA)
basic 14	DL-p-Chloroamphetamine HCI	0 resolved	3	7



Bringing selectivity with extended solvent range

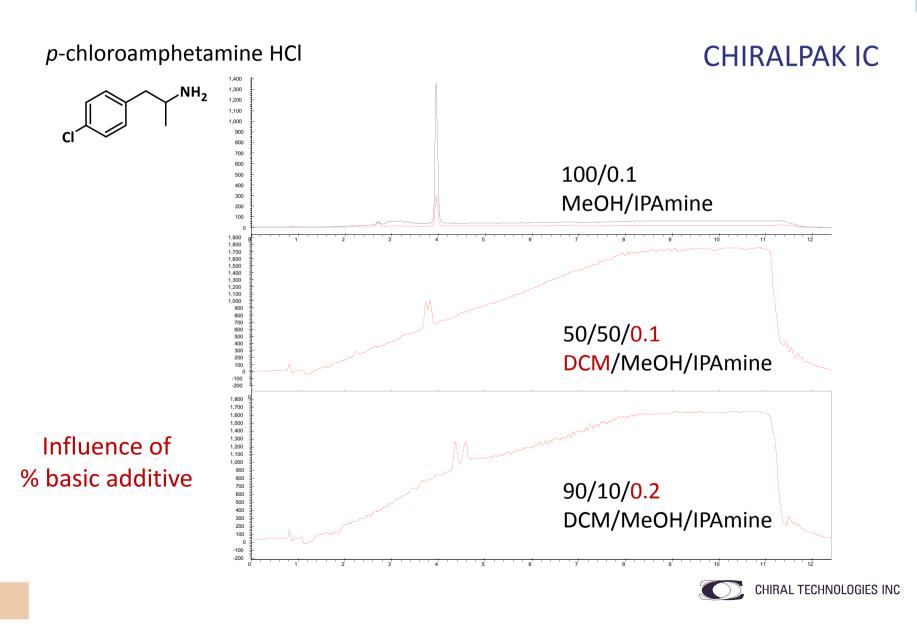


The success rate

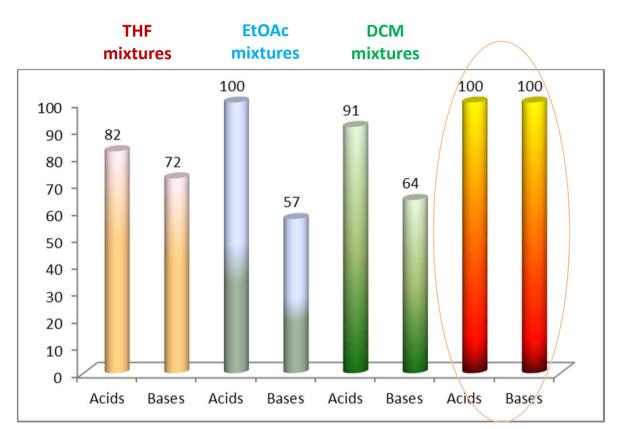

Global success extended solvent range

- ✓ All molecules resolved in the extended solvent range
- ✓ For acidic compounds, EtOAc leads to higher success rate
- ✓ For bases, it may be important to further investigate higher % basic additive

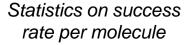




Overview of conditions tested - THF

THF / MeOH
90 /10
+
0.2 additive

- ✓ Proper elution strength (>70%)
- ✓ Resolution success rate:
 - ✓ 82% for acids
 - ✓ 72% for bases

Diperodon

CHIRALPAK IA THF/MeOH/IPAmine

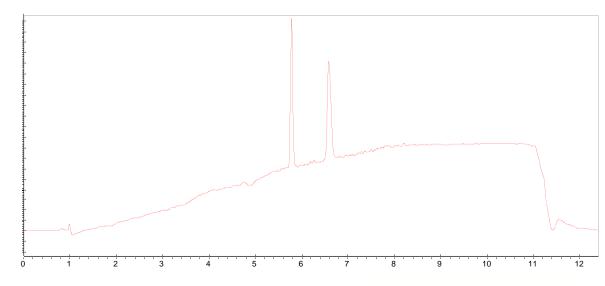




Overview of conditions tested - EtOAc

EtOAc / MeOH 90 /10 + 0.2 additive

- ✓ Proper elution strength (>70%)
- ✓ Resolution success rate:
 - ✓ 100% for acids
 - ✓ 57% for bases
- ✓ For bases, the addition of more alcohol and higher % IPAmine should be investigated

Verapamil

CHIRALPAK IB EtOAc/MeOH/IPAmine

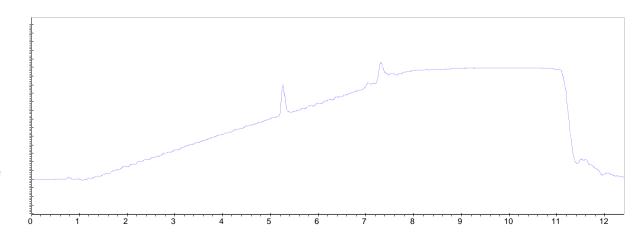




Overview of conditions tested - DCM


DCM/ MeOH 50 /50 or 90/10 + 0.2 additive

DCM mixtures

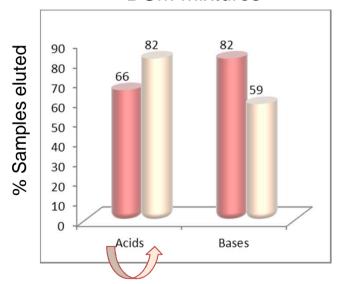

From 50% MeOH to 10% MeOH in DCM

- Different tendencies for acids and bases
 - ✓ Acids elute better and reach higher success rates with 10% alcohol
 - ✓ Bases need more alcohol for elution and resolution
- ✓ Resolution success rate:
 - ✓ 91% for acids
 - ✓ 64% for bases
- ✓ For bases, the addition of more alcohol and higher
 %IPAmine should be investigated

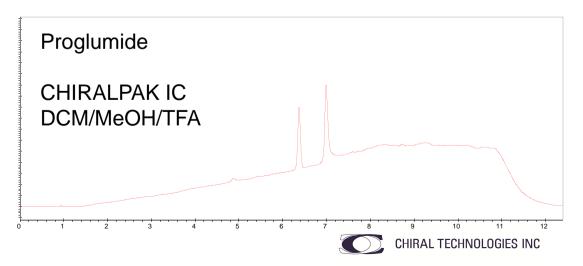

Column Performance Comparisons

Total 25 compounds (acids and bases)					
Co-solvents	IA	IB	IC	ID	ΙE
95% THF/5%MeOH	10	6	14	12	7
90% EtOAc/10%MeOH	13	11	13	12	9
90% DCM/10%MeOH	11	7	9	8	8
Total Hits	34	24	36	32	24
Co-solvents					
95% THF/5%MeOH					
90% EtOAc/10%MeOH IC = IA > ID > IB > IE					
90% DCM/10%MeOH		E > IB			
Totals IC > IA > ID > IE = IB					



Conclusions

- ✓ The use of extended solvent range in SFC for chiral applications offers advantages to:
 - ✓ Increase solubility of the samples
 - ✓ Avoid instability of molecules sensitive to alcohols
 - ✓ Broaden selectivity profiles and increase success rate

 High interest to explore all analytical options of SFC now that new powerful SFC equipment is available
- ✓ Screening strategies should be adapted accordingly in terms of co-solvent composition and presence of additives
- ✓ Other solvents (MtBE, ...) and conditions (other additive combinations) could be further explored



Acknowledgements

- Frank Riley, Tony Yan and other scientists at Pfizer (Groton, CT)
- ➤ Mark McDonald of Chiral Technologies Inc

